


Lymphosep 1077 (Histopaque)

Cat. N.	Product	Size	Price
IS-1077-500	Lymphosep 1077 (Histopaque)	1 x 500 ml	€ 95
IS-1077-600	Lymphosep 1077 (Histopaque)	6 x 500 ml	€ 450
IS-1077-100	Lymphosep 1077 (Histopaque)	6 x 100 ml	€ 168

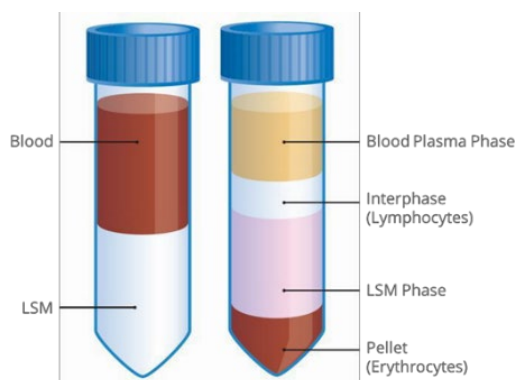

Suitable for:

- The simple, rapid isolation of lymphocytes from whole blood
- Peripheral blood
- Isolation of Mono Nuclear Cells (MNC)
- Whole blood, diluted that has been treated with anti-coagulants or defibrinating agents.
- Lymphosep is a Lymphocyte Separation Media / Histopaque

TABELLA COMPARATIVA

Immunological Sciences vs. Sigma-Aldrich

Immunological Sciences			Sigma Aldrich		
Cat.n.	Size	Price	Price	Cat.n.	Size
IS-1077-500	1 x 500 ml	€ 95	€ 199	10771-500ML	1 x 500ml
IS-1077-100	6 x 100 ml	€ 168	€ 185	10771-6X100ML	6 X 100 ML

CONDIZIONI DI VENDITA:

Prezzi: al netto dell'IVA (22%)

Consegna Pronta (salvo il venduto)

Trasporto gratuito per ordini superiori a € 200

Offerta valida fino al 15/12/2026

Codice Offerta: LYMPHO-2026



SOCIETÀ ITALIANA CHIMICI
Life Sciences

Lymphosep 1077 (Histopaque)

Theoretical pH: 7.0 ± 0.5

Osmolality: 300 mOsm/kg ± 20

Density: 1.077 ± 0.001

Colour: colourless, clear solution

Storage conditions: Room temperature, protected from light

Shelf life: 24 months

Sterility tests:

- bacteria in aerobic and anaerobic conditions
- fungi and yeast

Endotoxin: < 10 EU/ml

Composition: Displayed on website and also available on request

Recommended use:

- Respect storage conditions of the product
- Do not use the product after its expiry date
- Store product in an area protected from light (not necessary for saline solutions).
- Manipulate the product in aseptic conditions (e.g. : under laminar air flow)
- Wear clothes adapted to the manipulation of the product to avoid contamination (e.g. : gloves, mask, hygiene cap, overall...)

The product is intended to be used for in vitro diagnostic and scientific purposes. Do not use it in therapy, human or veterinary applications.

Applications:

Lymphosep 1077 is designed for the simple, rapid isolation of lymphocytes from whole blood , peripheral blood , to isolate Mono Nuclear Cells (MNC)

The product is suitable for whole blood , diluted that has been treated with anti-colagulants or defibrinating agents.

For best results, use blood drawn less than 2 hours before. Do not use blood more than 24 hours from when it was drawn.

Lymphosep 1077 is also called Histopaque 1077

PROTOCOL OF USE:

- 1) Thoroughly mix the Lymphosep 1077 by inverting the bottle gently.
- 2) Aseptically transfer 3 ml of Lymphosep 1077 to a 15 ml centrifuge tube.
- 3) Mix 2 ml of defibrinated or heparined blood with 2 ml of physiological saline (PBS w/o Ca w/o Mg) or balanced salt solution .
- 4) Carefully layer the diluted blood over 3 ml of Lymphosep (room temperature) in a 15 ml centrifuge, creating a sharp blood- Lymphosep 1077 interphase. DO NOT MIX! The quality of the separation is dependent upon a sharp interphase between the lymphocytes and the solution.
- 5) Centrifuge the tube at 400G at room temperature for 15 to 30 minutes. Centrifugation should sediment erythrocytes and polynuclear leukocytes and band mononuclear lymphocytes above the Lymphosep 1077.
- 6) Aspirate the top layer of clear plasma to within 2-3 mm above the lymphocyte layer.
- 7) Aspirate the lymphocyte layer plus about half of the Lymphosep 1077 layer below it and transfer it to a centrifuge tube. Add an equal volume of buffered balanced salt solution to the lymphocyte layer in the centrifuge tube and centrifuge for 10 minutes at room temperature (18°C to 25°C) at a speed sufficient to sediment the cells without damage i.e., 160-260 g. Washing the cells removes Lymphosep 1077 and reduces the percentage of platelets.
- 8) Wash the cells again with buffered balanced salt solution (ISL0615) and resuspend in the appropriate medium for your applications.